

Care and Use of the VP 772FC-2



Figure 1. The VP 772FC-2 is designed to separate magnetic beads in 1-liter bottles.

OPERATION



1. Your 1-liter bottle should slide into the VP 772FC-2 easily. Some 1 liter's bottles will vary in other diameter. If your bottle does not fit, do not force it in. Please contact us for technical service.



2. Insert the bottle into the device. Please cap over the bottle and twist to lock into place. This will secure the bottle in the device.
3. Allow the magnetic beads to separate according to your protocol.
- 4.



5. Uncap the bottle, grab the handle on the device and decant the supernatant.

PRODUCT MAINTENANCE

GENERAL PRODUCT CARE

When not in use, place device away from ferrous materials.

The deck and case of the Magnetic Tumble Stirrer are made of PVC. To clean the deck, wipe down with a cloth and mild detergent followed by a water wipe. **Do not immerse The Magnetic Bead Separation Device in liquid.**

For chemical compatibility of ABS please see the following link for more information:

<http://www.coleparmer.com/techinfo/chemcomp.asp>

Caution

The VP 772FC-2 uses a large powerful magnet. Please use caution when placing is near ferrous materials as it will pull any magnetic material to it.

WARRANTY

V&P Scientific, Inc. warrants this product to be free from defects in material and workmanship when used under normal laboratory conditions for one year. This warranty begins from the date of delivery by V&P Scientific.

In the event this product fails under normal laboratory conditions within the specified period of time because of a defect in material or workmanship, V&P Scientific will, at its option, repair or replace the product. Damage to the product caused by user negligence is not covered.

Please keep the special shipping carton in case the unit needs to be shipped back to V&P Scientific. Please contact V&P Scientific at the address below for return authorization and shipping instructions.

This warranty is made in lieu of other warranties expressed or implied including the warranties of merchantability and fitness for a particular purpose. V&P Scientific shall not be liable for loss or damages arising from the use of these products nor for consequential damages of any kind.

If assistance is required, contact V&P Scientific, Inc. at 858-455-0643 or sales@vp-sci.com.